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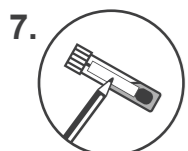
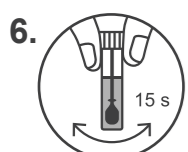
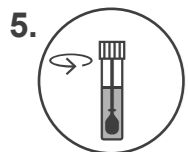
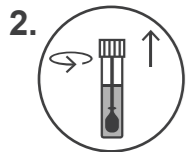
info@invitek.com
www.invitek.com

ENGLISH

Stool Collection Tube with DNA Stabilizer Instructions

RUO Research use only

Please read these instructions carefully before using the product.

INTENDED USE:

Collection of human stool samples for DNA stabilization, storage and extraction (PSP® Spin Stool DNA Plus Kit / PSP® Spin Stool DNA Basic Kit). The DNA is protected from degradation for up to 3 months at room temperature in the DNA Stabilizer. The product is intended for use under supervision of professionals only, such as technicians, physicians and biologists trained in molecular biological techniques.

CONTENTS:

Stool Collection Tube with integrated spoon in lid, label for donor description, each tube contains 8 ml of DNA Stabilizer reagent.

LIMITATIONS OF THE PROCEDURE:

Empty your bladder before beginning the collection. The sample size should NOT exceed the size of the spoon. Optimal is a sample size comparable to a small bean. When collecting liquid fecal sample transfer about 4 - 5 spoons of sample into the Stool Collection Tube.

IMPORTANT:

Before use make sure that the Stool Collection Tube has room temperature.

1. Collect fecal sample free of urine or toilet water. Toilet paper or tissues may be required.

2. Unscrew the lid of the Stool Collection Tube. Store the tube upright. Do NOT spill the stabilizing liquid in the tube.

3. Use the integrated spoon in the lid to collect a small stool sample. Do NOT collect too much sample.

4. Carefully transfer the spoon with the sample into the Stool Collection Tube with DNA Stabilizer.

5. Screw lid tightly onto the tube again.

6. Shake tube for 15 sec to mix fecal sample with DNA Stabilizer reagent.

7. Label the tube with donor name, birth date and collection date using provided label.

STORAGE:

Store at room temperature 15 – 30°C (59 – 86°F).

SHIPMENT:

Ship according to locally applicable guidelines and regulations such as UN 3373 (Biological Substance, Category B).

SAFETY PRECAUTIONS:

Do NOT use the Stool Collection Tube if liquid content of the tube is spilled.

Do NOT remove DNA Stabilizer reagent from the tube.
Do NOT ingest the contents of the tube.
Keep out of the reach of children. Wash with water if DNA Stabilizer reagent comes in contact with skin or eyes. Protect container from warming up and direct sunlight. Keep only in the original container.

WARNING:

Hazard statements

H319: Causes serious eye irritation.

H412: Harmful to aquatic life with long lasting effects.

Precautionary statements

P264: Wash hands, forearms and face thoroughly after handling.

P273: Avoid release to the environment.

P280: Wear protective gloves/protective clothing/eye protection/face protection/hearing protection.

P305+P351+P338: IF IN EYES: Rinse cautiously with water for several minutes. Remove contact lenses, if present and easy to do. Continue rinsing.

P337+P313: If eye irritation persists: Get medical advice/attention.

P501: Dispose of contents/container to hazardous or special waste collection point, in accordance with local, regional, national and/or international regulation.

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ALS Life Sciences Portugal, S.A.